

CONTROL III®

LABORATORY GERMICIDE

READY TO USE

**BACTERICIDAL
FUNGICIDAL
*VIRUCIDAL**

**REQUIRES NO ACTIVATORS
CONTENTS ARE READY FOR USE**

ACTIVE INGREDIENTS

n-alkyl
(60% C14, 30% C16, 5% C18, 5% C12)
dimethyl benzyl ammonium chloride
.....0.0781%

n-alkyl
(68% C12, 32% C14) dimethyl
ethyl benzyl ammonium chloride
.....0.0781%

INERT INGREDIENTS.....99.8438%

TOTAL..... 100.000%

**KEEP OUT OF REACH
OF CHILDREN**

CAUTION

SEE LEFT PANEL FOR ADDITIONAL
PRECAUTIONARY STATEMENTS

ONE GALLON (128 FL. OZ.)

EPA EST. #40873 CA-01/GA-01

EPAREG.#55364-4

MARIL PRODUCTS, INC.

15421 Red Hill Avenue, Suite D
Tustin, CA 92780 USA
800-546-7711 • 714-544-7711

DIRECTIONS FOR USE

It is a violation of Federal Law to use this product in a manner inconsistent with its labeling. This product is not to be used as a terminal sterilant/high level disinfectant on any surface or instrument that (1) is introduced directly into the human body, either into or in contact with the bloodstream or normally sterile areas of the body, or (2) contacts intact mucous membranes but which does not ordinarily penetrate the blood barrier or otherwise enter normally sterile areas of the body. This product may be used to pre-clean or decontaminate critical or semi-critical medical devices prior to sterilization or high-level disinfection.

DISINFECTION:

Control III Laboratory Germicide is a "Ready-To-Use" disinfectant which is designed for use as a disinfectant in hospitals, doctors' offices, dentists' offices, nursing homes, clinics and surgi-centers. It is formulated to disinfect hard, non-porous, inanimate surfaces such as counter tops, beds, instruments, fixtures and equipment.

1. Prior to disinfection with Control III Laboratory Germicide, all surfaces must be thoroughly cleaned to remove gross filth or heavy soil.
2. Use full strength. Do not dilute.
3. Apply Control III Laboratory Germicide with a clean sponge, cloth or by a trigger sprayer so as to wet all surfaces thoroughly. For disinfection, surfaces must remain wet for ten minutes. Allow to air dry.
4. Control III Laboratory Germicide

is classified as a low-level disinfectant.

OPERATING ROOM GERMICIDE:

Apply CONTROL III Laboratory Germicide solution to a clean sponge or cloth. Wipe down surgery tables, instrument stands, and other equipment so as to wet all surfaces thoroughly. Allow to air dry. (See Disinfection directions)

SURGICAL INSTRUMENT STORAGE:

Place instruments in clean storage tray container. Fill container with CONTROL III Laboratory Germicide solution to fully immerse instruments. Cover container. In normal use storage solution can be retained for thirty days. Replace solution more often if frequent instrument removals and additions are made. Instruments should be autoclaved after soaking in Control III prior to reuse. (See Disinfection directions)

Effective Against:

Pseudomonas aeruginosa (PRD-10); Staphylococcus aureus (ATCC 6538); Salmonella choleraesuis (ATCC 10708); Influenza A2 virus*; Herpes simplex type 1*; Adenovirus type #5; Vaccinia viruses*; Trichophyton mentagrophytes, on hard, non-porous, inanimate surfaces.

Note: Control III Laboratory Germicide has not been tested for effectiveness against Mycobacterium Tuberculosis and must not be relied upon when a Tuberculocidal product is desired.

HIV-1: See Technical Bulletin "Efficacy of Control III Products For the Control of the Human Immunodeficiency Virus Type 1 (AIDS VIRUS).

KILLS HIV-1 ON PRE-CLEANED ENVIRONMENTAL SURFACES/OBJECTS PREVIOUSLY SOILED WITH BLOOD/BODY FLUIDS in health care settings in which there is an expected likelihood of soiling of inanimate surfaces/objects with blood or body fluids can be associated with the potential for transmission of Human Immunodeficiency Virus Type 1 (HIV-1) (associated with AIDS).

SPECIAL INSTRUCTIONS FOR CLEANING AND DECONTAMINATION AGAINST HIV-1 OF SURFACES/OBJECTS SOILED WITH BLOOD/BODY FLUIDS. PERSONAL PROTECTION:

Clean-up should always be done wearing protective latex gloves, masks, and eye protection.

CLEANING PROCEDURES:
Blood and other body fluids

containing HIV must be thoroughly cleaned from surfaces and objects before application of this product.

CONTACT TIME:

Allow surface to remain wet for 10 minutes.

DISPOSAL OF INFECTIOUS MATERIALS:

Bloody body fluids, cleaning materials and clothing should be autoclaved and disposed of according to local regulations for infectious waste disposal.

PRECAUTIONARY STATEMENTS: HAZARDS TO HUMANS AND DOMESTIC ANIMALS**CAUTION:**

Causes moderate eye irritation. Avoid contact with eyes, skin, or clothing. Wash thoroughly with soap and water after handling.

STORAGE AND DISPOSAL:

Do not contaminate water, food, or feed by storage and disposal.

PESTICIDE STORAGE:

Store in a dry place no lower in temperature than 50°F or higher than 120°F.

PESTICIDE DISPOSAL:

Wastes resulting from the use of this product must be disposed of on site or at an approved waste disposal facility.

CONTAINER HANDLING:

Nonrefillable container. Do not reuse or refill this container. Offer for recycling if available or wrap container and put in trash.

FIRST AID

If in eyes: Hold eye open and rinse slowly and gently with water for 15-20 minutes. Remove contact lenses, if present, after the first five minutes, then continue rinsing eye. Call a poison control center or doctor for treatment advice.

If on skin or clothing: Take off contaminated clothing. Rinse skin immediately with plenty of water for 15-20 minutes. Call a poison control center or doctor for treatment advice.

If swallowed: Call a poison control center or doctor immediately for treatment advice. Have a person sip a glass of water if able to swallow. Do not induce vomiting unless told to do so by the poison control center or doctor. Do not give anything by mouth to an unconscious person.

If inhaled: Move person to fresh air. If person is not breathing, call 911 or an ambulance then give artificial respiration, preferably by mouth-to-mouth, if possible. Call a poison control center or doctor for further treatment advice.

Note to Physician: Probable mucosal damage may contraindicate the use of gastric lavage.

HOT LINE NUMBER

For emergency medical advice, call your local poison control center (1-800-222-1222) or doctor. Have the product container or label with you when seeking medical advice or treatment.

P/N 10005 Rev. A

{Emerging Viral Pathogens}

This product qualifies for emerging viral pathogen claims per the EPA's 'Guidance to Registrants: Process for Making Claims Against Emerging Viral Pathogens not on EPA-Registered Disinfectant Labels' when used in accordance with the appropriate use directions indicated below.

This product meets the criteria to make claims against certain emerging pathogens from the following viral categories:

- Enveloped viruses

<i>For an emerging viral pathogen that is a/an...</i>	<i>...follow the directions for use for the following organisms on the label:</i>
Enveloped virus	Adenovirus type 5

[Product name] has demonstrated effectiveness against viruses similar to [name of emerging enveloped virus] on hard, non-porous surfaces. Therefore, [product name] can be used against [name of emerging enveloped virus] when used in accordance with the directions for use against Adenovirus type 5 on hard, non-porous surfaces. Refer to the [CDC or OIE] website at [pathogen-specific website address] for additional information.

[Name of illness/outbreak] is caused by [name of emerging enveloped virus]. [Product name] kills similar viruses and therefore can be used against [name of emerging enveloped virus] when used in accordance with the directions for use against Adenovirus type 5 on hard non-porous surfaces. Refer to the [CDC or OIE] website at [website address] for additional information